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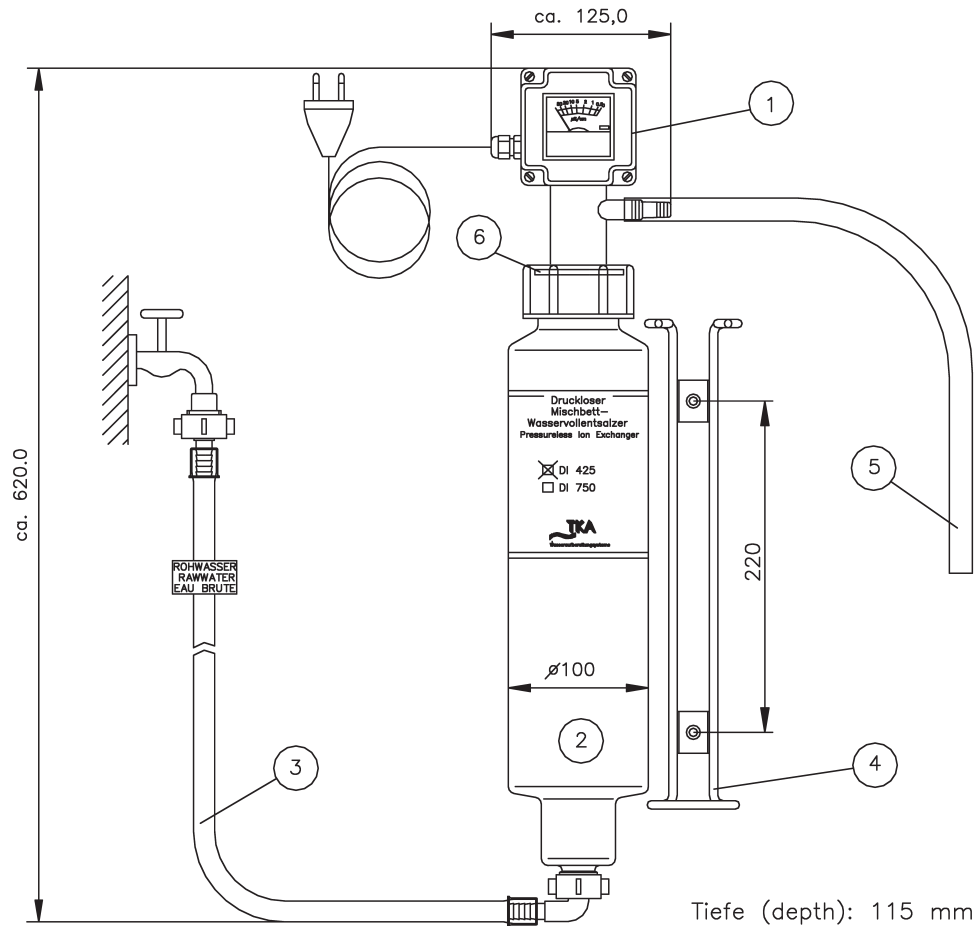
Art.-Nr.: 01.0425

Pressureless mixed bed ionexchange plastic cartridge

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Read these operating instructions carefully before installing and starting-up the system!

- | | |
|---|---|
| 1 Leitfähigkeitsmessgerät, analog für DI 425
Art.-Nr. 16.0001 | 1 Analog conductivity meter for DI 425
Art.-No. 16.0001 |
| 2 Drucklose Kunststoffpatrone
Art.-Nr. 01.0426 | 2 Plastic cartridge for pressureless operation
Art.-No. 01.0426 |
| 3 Rohwasserschlauch inkl. Dosierblende
Art.-Nr. 25.0036 | 3 Inlet hose incl. flow restrictor
Art.-No. 25.0036 |
| 4 Haltebügel
Art.-Nr. 18.0002 | 4 Bracket
Art.-No. 18.0002 |
| 5 Reinwasserschlauch ∅ 10 x 1,5 : 0,5 m
Art.-Nr. 18.0003 | 5 Outlet hose ∅ 10 x 1,5 : 0,5 m
Art.-No. 18.0003 |
| 6 Dichtung, ∅ 65 mm, weiß
Art.-Nr. 21.5013 | 6 Gasket, ∅ 65 mm, white
Art.-No. 21.5013 |

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Safety Precautions:

01. Using the cartridge in any other way than that described in these operating instructions invalidates the guarantee.
02. **Deionized water must on no account be consumed in any way. It is not fit for drinking!**

Mounting Instructions

01. Remove cartridge from bracket. Mount bracket near the point of use by using the enclosed screws and dowels.
02. Replace cartridge in bracket.
03. Connect the inlet hose with the **straight connector (flow restrictor integrated) to the water tap (connector is labelled "raw-water") and the angled connector to the unit.** Take care that the hose is not buckled.
04. Slide the outlet hose over the small outlet tube about half way up the conductivity meter (if necessary, warm the hose slightly). Connect the unit to the 230V outlet.
05. After opening of the tap, fully deionized water is supplied through the outlet hose.

ATTENTION: Never close off the hose connected to the purified outlet, or in any way impede free flow through it, as the plastic cartridge could burst under build-up pressure.

06. The unit can be used until the pointer of the conductivity meter leaves the green section or shows 20 $\mu\text{S}/\text{cm}$.

Change of Cartridge:

The cartridge is exhausted as soon as the pointer of the conductivity meter leaves the green marking or indicates 20 $\mu\text{S}/\text{cm}$.

01. Close water tap and remove electric plug from socket. Disconnect inlet hose from the cartridge. Let water run out of hose and the cartridge into water basin or a bowl until the cartridge is empty.
02. Remove cartridge spent from the bracket.
03. Unscrew the conductivity meter and dispose of the spent cartridge.
04. Fasten the inlet hose to the replacement cartridge and screw on the conductivity meter.
05. Fit the cartridge with the inlet hose in the bracket and fasten it.
06. Plug in the power plug. The unit is now ready for use.

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ECC-Declaration of conformity

According to EG Directive 89/336/EEC
- Electromagnetic Compatibility -

We hereby declare that the machine specified below conforms, in its design and construction, and in the versions we have introduced into the market, to the fundamental safety and health requirements of the EEC Directive 89/336/EEC.

This declaration loses its validity when changes which were not agreed to by us are made to the machine.

Designation of the machine: **Mixed-Bed Deionizer System for Purified Water**
Machine type: **TKA DI 425**
Article number: **01.0425**
Applicable EEC Directives: **EEC Directive Low voltage (73/23/EEC)**
EEC Directive Electromagnetic Compatibility (89/336/EEC)
Applied Standards: **EN 55014**

Niederelbert, 10 March 2003

Manufacturer / authorized representative

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Wall unit for direct connection to a water supply, with free flow through to the open purified water outlet. Supplies deionized water for laboratories and for doctors and dentists practices.

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The benefits:

- Supplies purified water at a very favourable low cost
- Easy to use
- Extremely effective
- Lightproof blue cartridge avoids growth of algae
- Wall mounting allows the purified water supply to be most practically positioned at the workplace

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Dimension, Performance and Specifications of TKA DI 425:

Flow rate, l/h	50
Capacity at 10°D TDS*, litres	425
Purified water quality, µS/cm	0,1 - 20
Max. feedwater temp., °C	30
Electrical requirements, V/Hz	230/50–60
Height** x diameter, mm	600 x 100
Approx. depth, mm	115
Approx. weight, kg	3
Order number	01.0425

* Total dissolved solids in the feedwater (10° D = 12.5°e)
** including conductivity meter

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Equipment supplied under the order number 01.0425

- Analog conductivity meter
- Cartridge filled with expendable resin
- Feed water hose incl. flow restrictor
(If you need to replace the feed water hose you are obligated to order the original TKA hose with the flow restrictor.)
- Purified water hose
- Wall mounting

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Carl Stuart Limited
TKA Wasseraufbereitungssysteme GmbH